


Operating instruction TUST37C and TUST38C



Simply insert the TUST37C (or TUST38C) into the phototube of your microscope, respectively with some LEICA microscopes in the **LeicaC1XTH port**. Next, fit the entire unit (LM digital adapter and camera) into the TUST37C (or TUST38C).

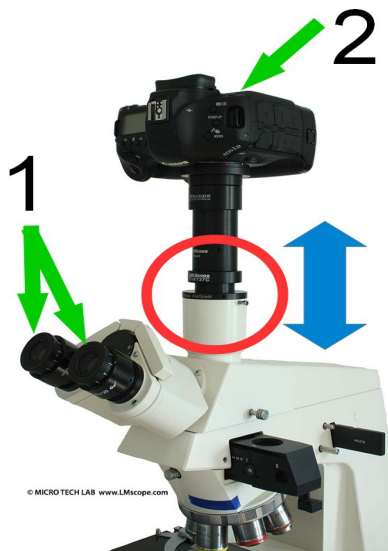

Then you should adjust parfocality (sharp image on observation tube and photo at the same time).

1) First, **focus the image** when looking through the **eyepicetubes**.

2) Now check, if the image is also in focus on the **switched-on camera (in Live View mode)**. If it's not, **loosen the knurled screw** (tube fixation) and **lift or lower** the camera unit until the image is as well sharp on the camera's display or the monitor.



If the optimal position has been found, please **fix the tube again** with the knurled screw.