



Operating Instruction: C-Mount Port 1x (TUST42C_SZ) for Olympus SZ Stereo - Microscopes

Set screw

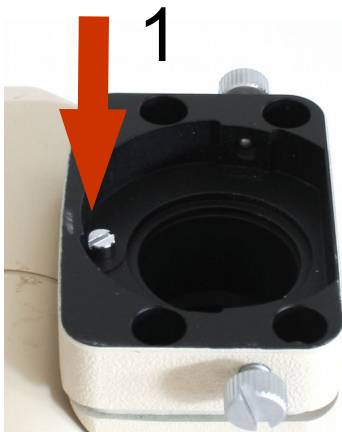

Remove the old adapter solution. Only the photo tube should remain. There is a small set screw on the photo tube, which is not needed with the new LM adapter solution. If the TUST42C_SZ can't be fitted properly because of this set screw, we recommend to simply remove the screw.

2

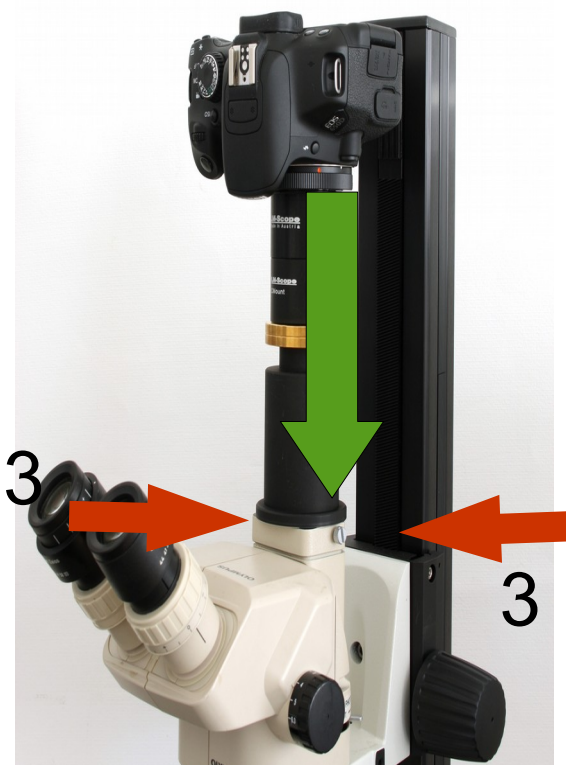

Remove the set screw (image 1). Now loosen the two knurled screws (3) and plug the TUST42C_SZ in to the photo tube (2). Then fix the C-Mount Port 1x by tightening the two knurled screws. The TUST42C_SZ should be mounted securely.