

Operating Instruction Nikon VT tube



After completing the installation of the entire camera unit (camera and LM digital adapter), an optical comparison between eyepicetube und phototube should be performed. Only by adjusting the parfocality (simultaneously focused image in the observation tube and on the photo) an ergonomic working is possible.

First the image is focused by looking through the eyepicetubes.

Now check, if the image is also well focused in the Live View mode of the camera (switched on). If this is not the case, the knurled screw (tube's fixation) has to be loosened and the camera unit has to be lifted or lowered until the image is also in focus on the camera's display or a monitor.

After finding the optimal location, the tube has to be fixed again with the help of the knurled screw.