Micro Tech Lab

www.LMscope.com

Installation instruction for motorised Stackshot Macro Rails by Cognisys on a LM stand

By means of the LM Mounting plate with M8-thread the Stackshot Macro Rail can be fixed easily on the LM Stand.




First of all remove the LM objective and the whole lens mounting unit from the LM stand.



After removal you will see the position pins on the stand rail.



Now place the LM assembly plate on the stand rail. Please take care that the LMScope label is at the bottom and the position pins are clicked into place.

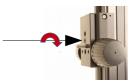

Fix the plate with the hex screw.



Place the Stackshot Macro Rail centrally on the M8setscrew and fix the unit with the washer and the M8-nut.



Check the axial alignment of the Stackshot Macro Rail with a small spirit level, then tighten the M8-nut.



Please check if the mounted Macro Rail can pass the screw smoothly before start-up.



Finally mount the camera and the objective unit again.