

Operating Instruction for tube adapters with 30 mm external diameter for LM digital adapter

	<p>Depending on the brand and model of the microscope, the interior diameter of the eyepiece tube might differ by a few tenths of a millimetre from the measurement required. For this reason, the tube adapter [TUST30W and TUST30S] comes with several different O-rings included. This allows the user to achieve an exact fit.</p>
	<p>Check how much extra space there is around the tube adapter without O-rings, and then use the larger or the smaller rings accordingly. You can also use a combination of larger and smaller O-rings.</p>
	<p>The tube adapter should fit into the eyepiece tube without being forced, and should not wobble when inserted.</p>
	<p>Next, fit the LM digital SLR universal adapter for 23.2 mm tubes into the TUST30 tube adapter and secure it in place by tightening the three Allen head screws.</p>
	<p>To achieve optimum focus, you have the possibility (as described above) of adjusting the distance between tube adapter and LM digital adapter using an Allen key.</p>